

Driver Installation

(For all Tecella products, except for the Pico)

Document Revision 1.1
December 23, 2020

Tecella
97 Chaumont Cir
Foothill Ranch, CA 92610, USA
www.tecella.com
info@tecella.com
+1-714-641-1709
+1-714-641-1569 Fax


1. Download and run the driver installer

For 64-bit systems:


http://downloads.tecella.com/driver/Driver_64bit_4.5.6.exe

For 32-bit systems:


http://downloads.tecella.com/driver/Driver_32bit_4.5.6.exe

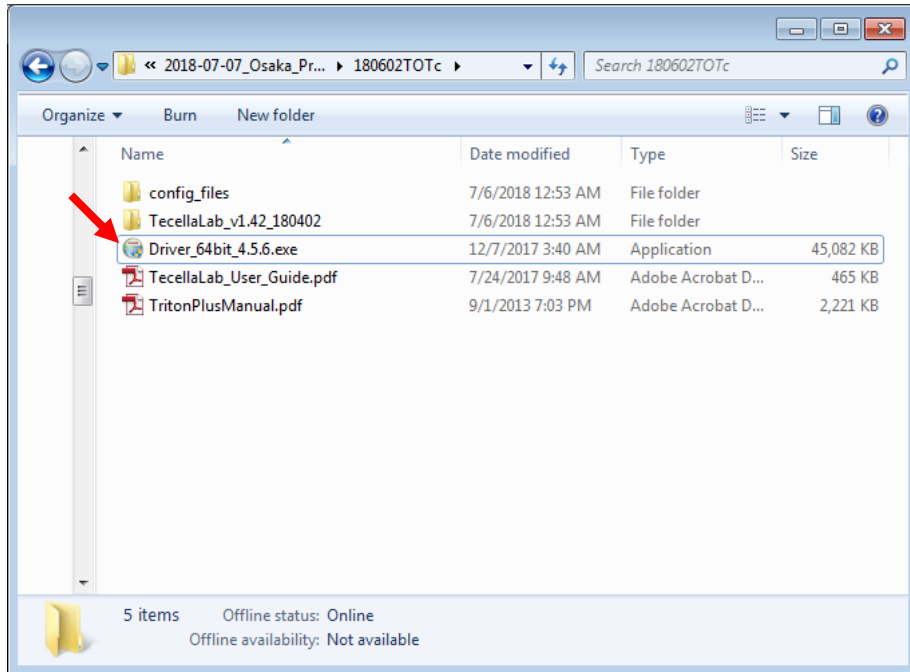


2. Click on “Next”

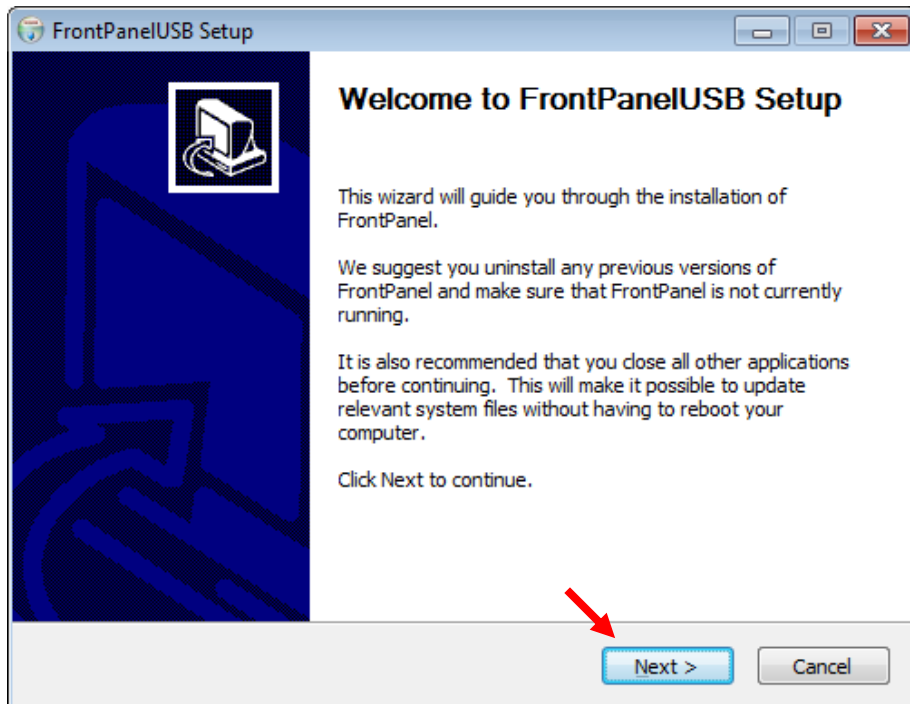


3. Click on "I Agree"

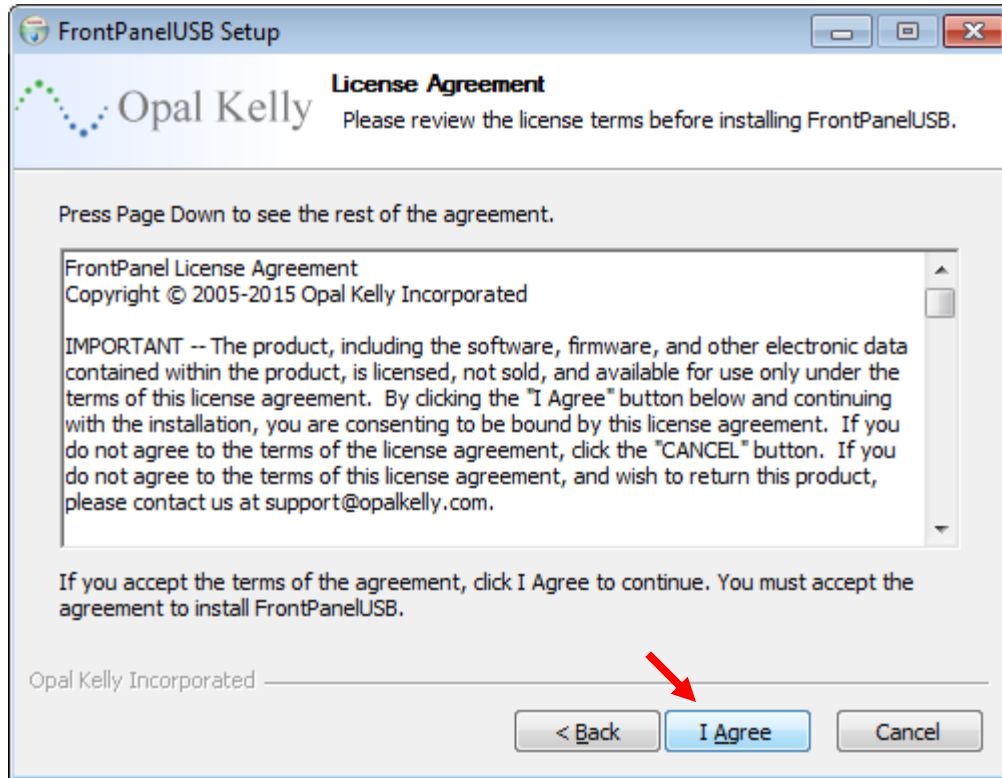

4. Uncheck everything except "FrontPanel USB Driver".
"FrontPanel USB Driver" is checked by default, and cannot be unchecked.

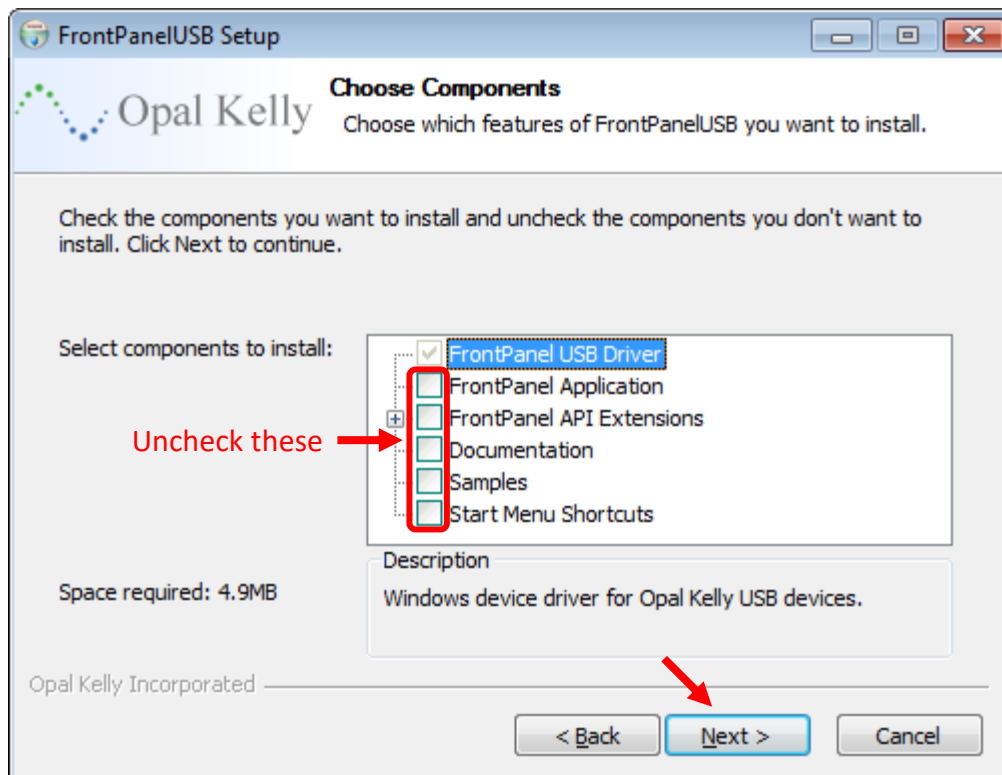

5. Click on "Install"

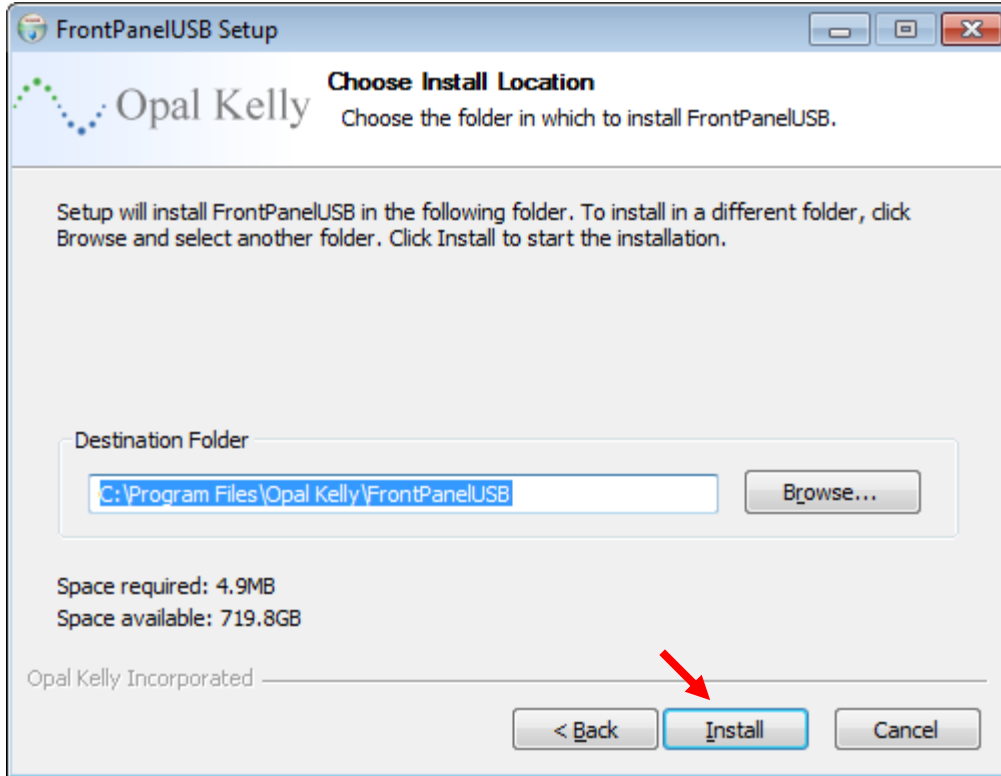

6. Click on "Finish" to complete driver installation

